

TSK-GEL® CM-3SW Products

| | | |
|----------------------------|-------------------------------------|---------------------|
| Part Numbers: | 07162, 7.5mm ID X 7.5CM, 10 μ m | 07650, Guardgel Kit |
| Counter Ion: | Na ⁺ | |
| Small Ion Capacity: | > 0.3 meq/gram | |

This sheet contains the recommended operating conditions and the specifications for TSK-GEL CM-3SW columns. Installation instructions and column care information are described in a separate Instruction Manual.

A. OPERATING CONDITIONS

1. Shipping Solvent: Methanol (Replace the shipping solvent with water before changing to the mobile phase of choice.)
2. Max. Flow Rate: 1.2 mL/min
When a buffer with high viscosity is used, the maximum flow rate may have to be reduced so as not to exceed the maximum pressure drop. When changing solvents, use a flow rate equal to 50% of the maximum flow rate.
3. Standard Flow Rate: 0.5 - 1.0 mL/min
4. Max. Pressure: 20 kg/cm² = 375 psi
5. pH Range: 2.0 - 7.5
6. Salt Conc.: \leq 0.5 M
7. Organic Conc.: 0 - 100% for aqueous soluble organic solvents. Make gradual solvent changes using a shallow gradient at low flow rate.
8. Temperature: 10 - 45°C. Reduce flow rate when operating below 10°C.
9. Cleaning Solvents: (1) Buffer with 0.5 - 1.0M salt, or
(2) pH 2 - 3 buffer, or
(3) Buffer containing methanol or acetonitrile, or
(4) Urea or nonionic surfactant in buffer.
NOTE: Choose a cleaning buffer based on sample properties, e.g. use (1) to remove strongly adsorbed proteins, (2) to reduce the negative charge on the biopolymer, (3) to decrease hydrophobic adsorption, and (4) to increase protein solubility.
10. Storage: Store the column in methanol at ambient temperature when it will not be used the next day. For overnight storage flush the column with the low salt concentration mobile phase at 0.2 mL/min. Prevent air from entering the column!
11. Column Protection: The use of guard columns is recommended to prolong the life of the analytical column. Guard column life depends greatly on sample cleanliness. As a general rule, guard columns should be replaced after every 30-40 sample injections, when the peaks become excessively wide, or when the peaks show splitting.

B. SPECIFICATIONS

The performance of TSK-GEL CM-3SW columns is tested under the conditions described in the Data Sheet. All columns have passed the following quality control specifications:

1. Number of Theoretical Plates (N): \geq 1,300
2. Asymmetry Factor (AF): 0.8 - 1.6

DS1032 Revised 24JANUARY2003

Tosoh Bioscience LLC
3604 Horizon Drive, Suite 100
King of Prussia, PA 19406
Phone: (484) 805-1219
Orders and Technical Service: (800) 366-4875
FAX: (610) 272-3028

Tosoh Bioscience GmbH
Zettachring 6
D-70567 Stuttgart
Germany
Phone: 0711/13257 0
FAX: 0711/13257 89